




KONICA MINOLTA

Spectrophotometer CM-3630
PaperControl Software

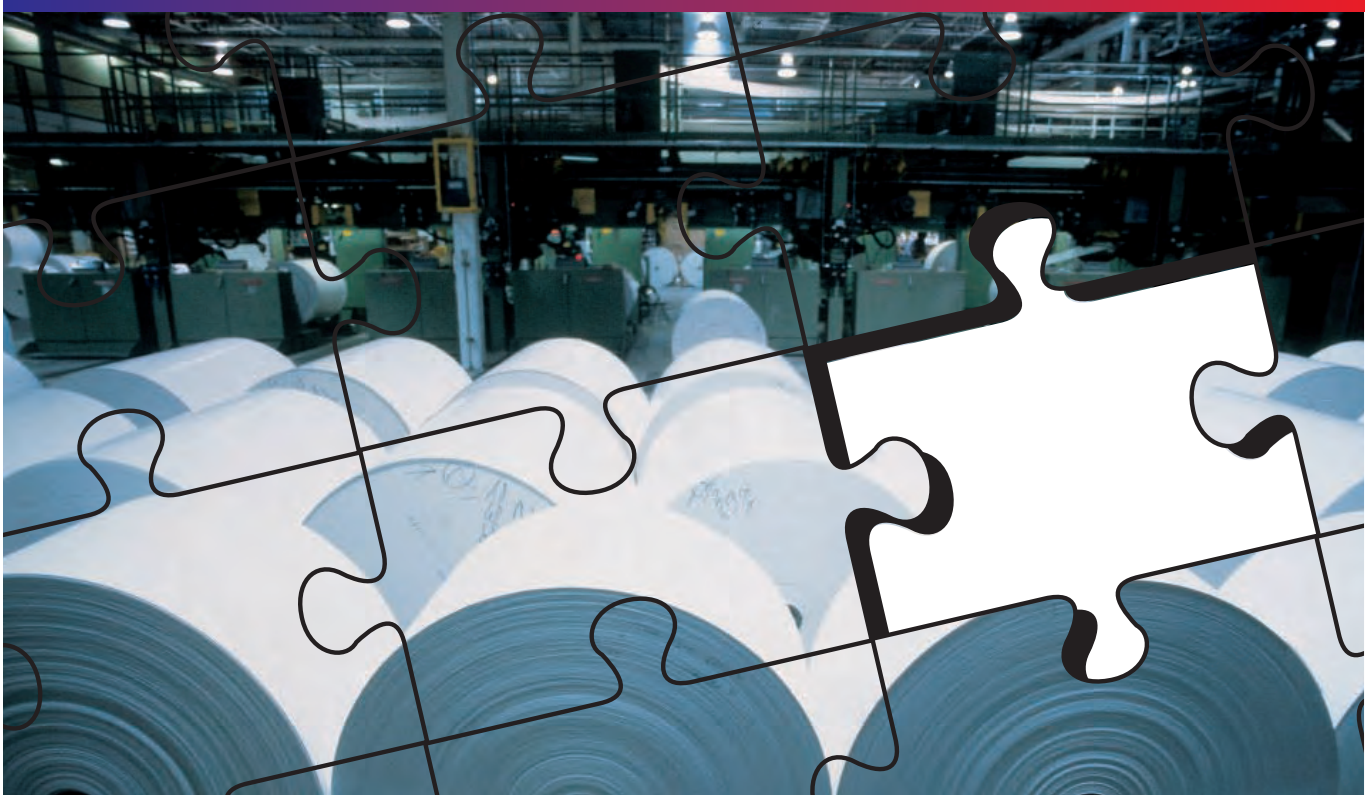

The missing link to total quality
Solutions made to measure the
paper industries' needs.

Introducing Konica Minolta's high-efficiency ISO
solution
for precise whiteness and color measurement

The essentials of imaging




Exactly what you've
been waiting for:
Introducing our state-of-the-
art Spectrophotometer
CM-3630 and PaperControl
software for the edge in
quality control.



Highly competitive markets, such as the paper, pulp and cellulose industry, require a commitment to top quality throughout. With Konica Minolta's state-of-the-art Spectrophotometer CM-3630 and the accompanying PaperControl software you can precisely monitor production at all times.

Designed to keep an eye on the optical properties of pulp and paper during the complete production process, the CM-3630 measures brightness, opacity, fluorescence, color, whiteness and tint. To meet your exacting standards, PaperControl software has been developed in close cooperation with the paper, pulp and cellulose industry.

Finally a professional solution to ensure uncompromised quality

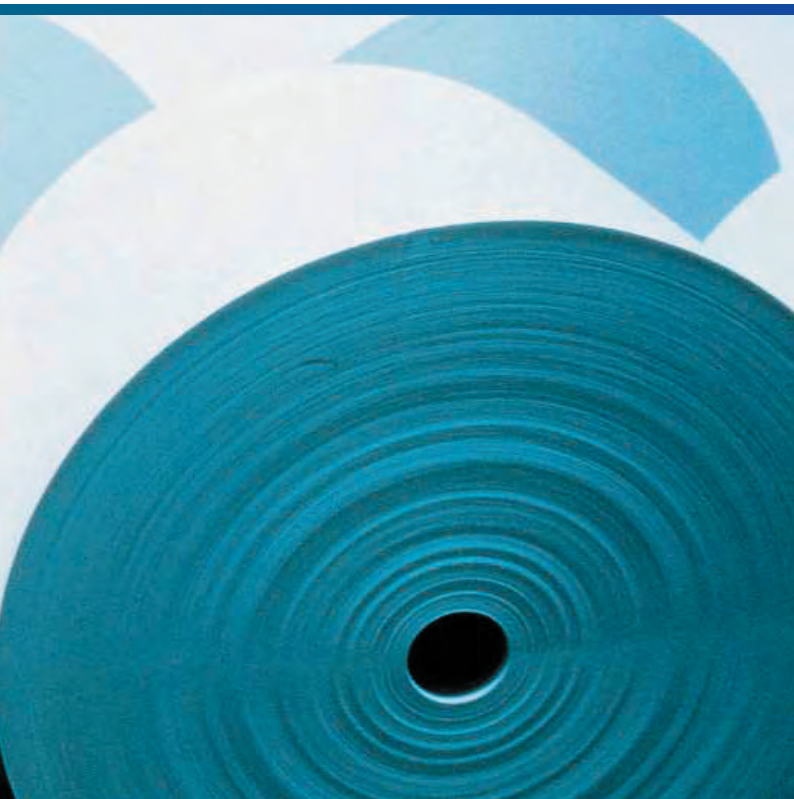

Your professional quality agent right on the production line: Uncompromised quality thanks to advanced PaperControl software and perfect adherence to ISO standards.

Measuring Color Difference

In addition to production control and whiteness measurement, PaperControl also provides a Color Difference module, which displays the measurement readings as numerical data, spectral reflection curves or color space plots.

Quick Opacity Measurement

Measuring opacity can be done in a matter of seconds: Simply position the opacity jig on the white side and measure. Then rotate the opacity jig to the black side, measure again – that's it!

Patented* Numerical UV Control

Numerical UV Control (NUVC) makes calibration and adjustments to whiteness and tint a swift and quick procedure. In fact, the whole process is much faster than current methods used in conventional instruments. Since it's all done digitally, exact and reliable results are assured. And with the total absence of moving parts in the spectrophotometer's optical system, it is absolutely robust and completely maintenance-free! With patented NUVC, Konica Minolta has turned the once time-consuming and error-prone calibration process into a quick, accurate and reliable procedure.

Data compatibility

The CM-3630's numerical UV control system combined with its user calibration option allows easy adjustment of the instrument parameters to your current measurement system. Such precision guarantees that your brightness measurements will perfectly correspond with the data you've collected over the years.

*US Patent No 5,636,015.

ISO standards

The CM-3630 ensures adherence to ISO optical standards and a wide array of national standards worldwide. It provides exact conformity with ISO standards 2469 and 2470 for measuring ISO brightness. With PaperControl software, brightness, opacity, whiteness and yellowness indices are calculated and displayed in accordance with several international and national standards such as: ISO, SCAN, DIN, TAPPI, CPPA, and AFNOR.

Whiteness Measurements

To get correct measurement readings of paper with Optical Brightening Agents, the amount of UV radiation must be controlled and adjusted. PaperControl offers several methods to adjust the UV radiation. To prevent the triplet effect exhibited by some special Optical Brightening Agents, the CM-3630 allows you to switch the light source between full power and soft flash mode.





Used by STFI
as reference instrument
for measuring white paper standards

Konica Minolta - a synonym for reliability and expertise throughout the paper industry


Konica Minolta manufactures its advanced measuring instruments with meticulous care. Strict quality controls ensure that each product rigorously conforms with exacting industry standards. For over 60 years, Konica Minolta has offered innovative, affordable solutions that deliver solid value through their reliability, durability, accuracy, precision and repeatable performance.

Technical Data

Model	CM-3630
Illumination/viewing system	Reflectance:d/0 (diffused illumination, 0-degree viewing) Conforms to ISO 2469, JIS P8148, DIN 53145-1 and DIN 53145-2 standards.
Light-receiving element	Silicon photodiode array (dual 40 elements)
Spectral separation device	Diffraction grating
Wavelength range	360nm to 740nm
Wavelength pitch	10nm
Reflectance range	0 to 200%; resolution: 0.01%
Light source	Pulsed xenon lamps (x 3)
Measurement time	Approx. 1.5 seconds (for measurements of fluorescent colors, at 9600 bps)
Minimum interval between measurements	Approx. 4 seconds; when reflectance measured, approx. 5 seconds; when fluorescent color
Measurement/illumination area	30mm/34mm
Inter instrument agreement	Mean ΔE^*ab 0.2 based on 12BCRA Series II color titels compared to values measured with master body
Repeatability	Spectral reflectance: Standard deviation within 0.1% Colorimetric values: Standard deviation within ΔE^*ab 0.02 (condition; white calibration plate measured 30 times at 10-second intervals)
Temperature drift	Spectral reflectance: Within +/- 0.10%/°C Color difference: Within ΔE^*ab 0.05/°C
UV adjustment	Instantaneous numerical adjustment
Control method	Directly connected to a computer
Interface	RS-232C format
Power	100-240V AC, 50-60Hz 25W AC (with a dedicated AC adapter)
Size (WxHxD)	300 x 585 x 315 mm
Weight	15.5 Kg
Operating temperature/humidity range	13 to 33°C, relative humidity 80% or less (at 30°C) with no condensation * Operating temperature/humidity range of products for North America : 13 to 33°C, relative humidity 80% or less (at 31°C) with no condensation
Standard accessories	White calibration plate CM-A 133, Zero calibration box CM-A119, AC Adapter AC-A12, AC Adapter AC-A21 (for Korea), RS-232C cable IF-A12, Accessory Case CM-A117, Dust Cover CM-A118, Unit Driver CM-A108
Optional accessories	Color Data Software PaperControl for CM-3630, Opacity jig CM-A134, RS-232C cable IF-A10, IF-A11, IF-A13, IF-A14, IF-A15
Specifications subject to change without notice	



Peace of mind for the pulp
and paper industry
Swift performance, reliable data
- everything under control

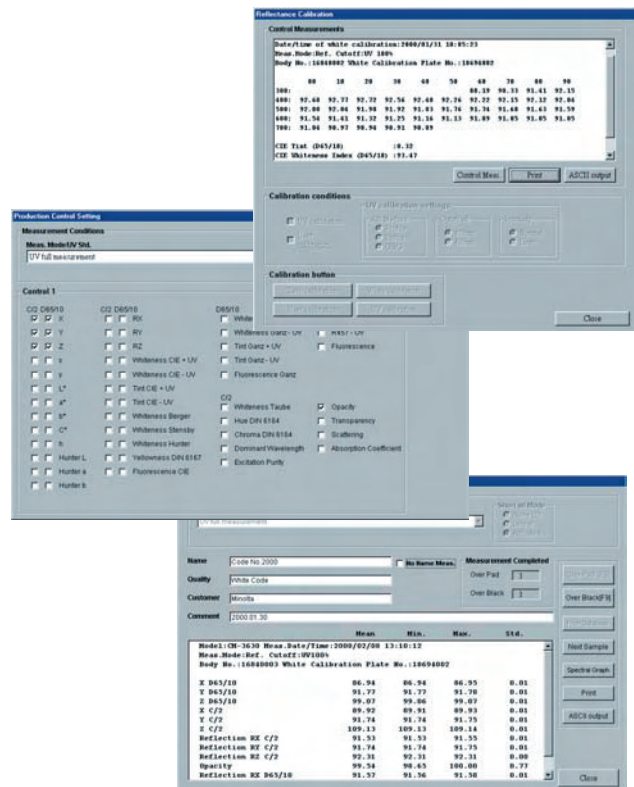

Production Control Module

PaperControl makes routine production control much easier, faster, more comfortable and reliable - in short: more professional. No more confusion by mixing production control and general QC tests as in conventional multi-purpose software. PaperControl is all you need to control your production line. All possible measurement parameter settings are already preprogrammed – for quick selection by the supervisor.


The production control module allows programming and labelling for a large number of jobs. Up to 9 different processes can be displayed in easy-to-understand menus. Thus, the operator gets valuable assistance in performing the most commonly used measurement routines. The end result is overall increased productivity.

The measurement data can be transferred to the production control process system. This ensures that specific test runs are being done under the same conditions – providing consistent, reliable results. And to make it even easier, all settings and measurements of this Windows-based software (Windows®95, Windows®98, and Windows NT®4.0) are keyboard-controlled.

Automatic calibration

Calibration is performed automatically for every conceivable combination of measurements in a few seconds. This valuable feature saves time and ensures reliable measurement data when new routines or tests are selected.

Windows® is a registered trademark of Microsoft Corporation, USA.

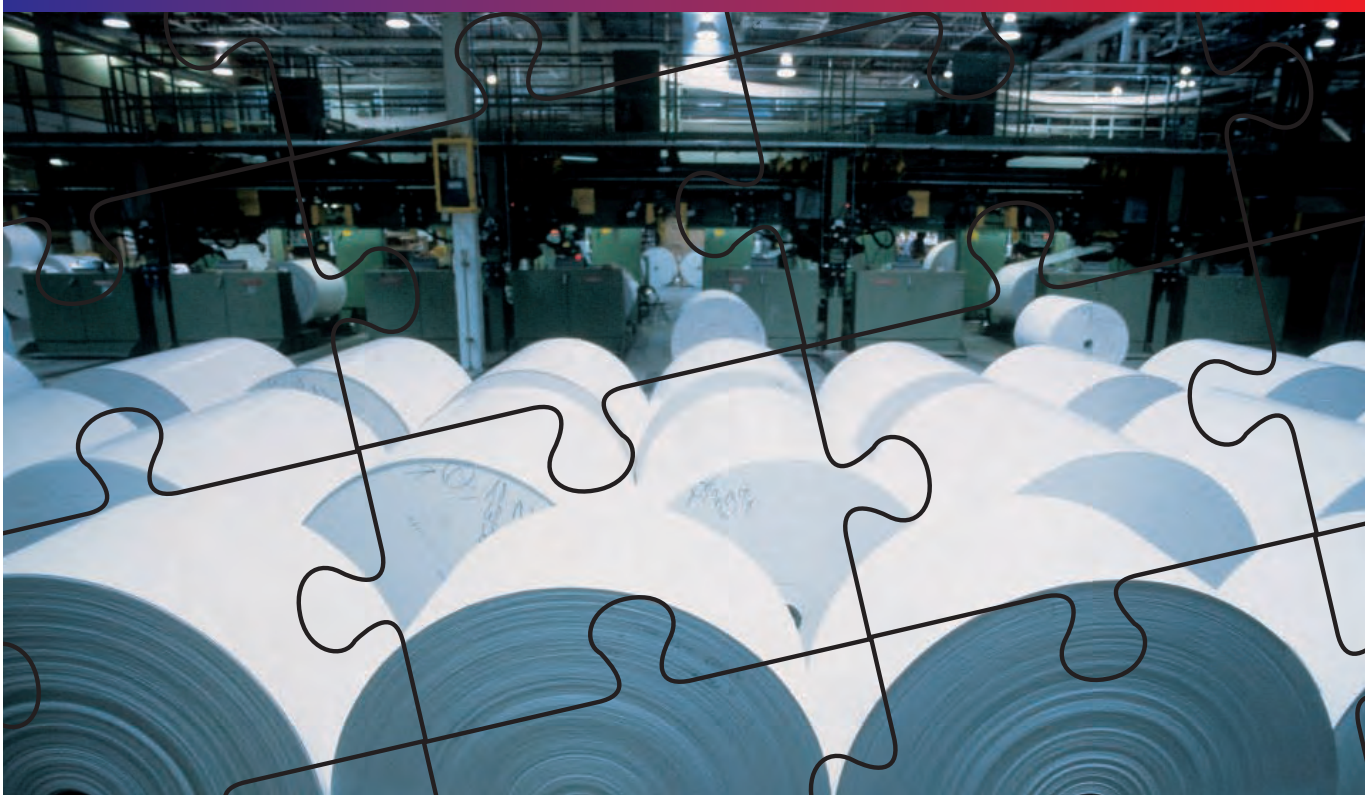

SAFETY PRECAUTIONS

For correct use and for your safety, be sure to read the instruction manual before using the instrument.



- Always connect the instrument to the specified power supply voltage. Improper connection may cause a fire or electric shock.



Certificate No : YKA 0937154
Registration Date : March 3, 1995



Certificate No : JQA-E-80027
Registration Date : March 12, 1997

KONICA MINOLTA SENSING, INC.
Konica Minolta Sensing Americas, Inc
Konica Minolta Sensing Europe B.V.

Osaka, Japan
New Jersey, U.S.A.
European Headquarter /BENELUX
German Office (International)
German Office (Germany)
French Office
UK Office
Italian Office
Swiss Office
Nordic Office
Austrian Office
Polish Office

Konica Minolta (CHINA) Investment Ltd.
SE Sales Division
SE Beijing Office
SE Guangzhou Office

Konica Minolta Sensing Singapore Pte Ltd.

KONICA MINOLTA SENSING, INC. Seoul Office

SE Sales Division	Shanghai, China	Phone : +86-021-5489 0202	Fax : +86-021-5489 0005
SE Beijing Office	Beijing, China	Phone : +86-010-8522 1551	Fax : +86-010-8522 1241
SE Guangzhou Office	Guangzhou, China	Phone : +86-020-3826 4220	Fax : +86-020-3826 4223
	Singapore	Phone : +65 6563-5533	Fax : +65 6560-9721
	Seoul, Korea	Phone : 02-523-9726	Fax : 02-523-9729

Addresses and telephone/fax numbers are subject to change without notice. For the latest contact information, please refer to the KONICA MINOLTA SENSING Worldwide Offices web page (link below).

<http://konicaminolta.com/about/se/contact.html>